

Viewing Images with Zeiss LSM Image Browser

1. Download folder and unzip as per Using PHPfileNavigator protocol. Make sure database *.mdb files are downloaded with the image files.
2. Open Zeiss LSM Image Browser.
3. Click **Open**.
4. Double click file you want to open.
5. Double click the database .mdb file you wish to open.
6. When window opens click the **Gallery** button.
7. Select Image you want and double click to open it.
8. Select how you want the image to appear: overlay (xy) or each channel plus overlay (split xy).
9. Click **Export** Button on Zeiss LSM Image Browser Menu
10. For Image type. If your image is:
xy: Select full resolution image - series
split xy: Select contend of image window - series
11. Save image as a **Tagged Image File [* .tif]**. Note: if you choose a 12 or 16 bit .tif file option you may have trouble opening it up depending on the program you choose to open it up in. It's best to save it as just a .tif.
12. Double Click .tif file to open it.

Images can now be opened in Photoshop or put into PowerPoint presentations.